

Microscope image measurement with QColor Camera

Standard Operating Procedure

Faculty Supervisor: Prof. Robert White, Mechanical Engineering (x72210)

Safety Office: Peter Nowak x73246 (Just dial this directly on any campus phone.)

(617)627-3246 (From off-campus or from a cell phone)

Tufts Emergency Medical Services are at x66911.

For more information on the program ImageJ and on plug-ins for the software see:

<http://rsb.info.nih.gov/ij/>

Revised: October 16, 2007

1.0 Material Requirements:

- 1.1 Equipment:** Fisher brand microscope with attached color CCD camera, Dell computer with video capture card and "Image J" software
- 1.2 Personal Protective Equipment:** Nitrile gloves and eye shields (standard cleanroom protective equipment to be worn at all times while in the cleanroom)

2.0 Procedure:

Take image:

1. Take an image from microscope as earlier described in the SOP "Microscope Image capture".
2. Save image (recommendation: save the image with the magnification factor in the filename (e.g. 5x_great_MEMS_device_1.jpg).

Measure distance:

1. Start ImageJ (microscope icon on desktop).
2. Open your previously saved image (File → Open...).
3. In order to set the pixels /distance ratio, select "Analyze" → "set scale". Enter "distance in pixels" and "know distance" according to the objective used to take the picture (5x, 10x, 20x or 40x). The conversion ratios are found in Table 1 below. Turn "Global" checkbox on and set units to um.
4. In order to measure, left click on the "Straight line selections" button in the toolbar.
5. Locate cursor to the starting point of measurement → left click, hold and drag to the desired end point of measurement. (Note: hold down shift to force the line to be at 90 degrees to the screen)
6. Select "Analyze" → select "Measure" → window with quantities pops up. (shortcut: Ctrl-m)

Comments:

1. ImageJ offers a number of other functions (e.g. label...), it is encouraged to further explore the program. Try "Analyze/Tools/Scale bar" to add a scale bar to your image.
2. ImageJ offers a lot of plug-ins for a variety of image analysis problems (e.g. contact angle measurement...), please visit the website.

Objective	Distance in Pixels	Known Distance (um)	Aspect Ratio
5x	0.822	1	1
10x	1.59	1	1
20x	3.20	1	1
40x	6.24	1	1

These parameters are for the QColor CCD camera with 1x1 binning.